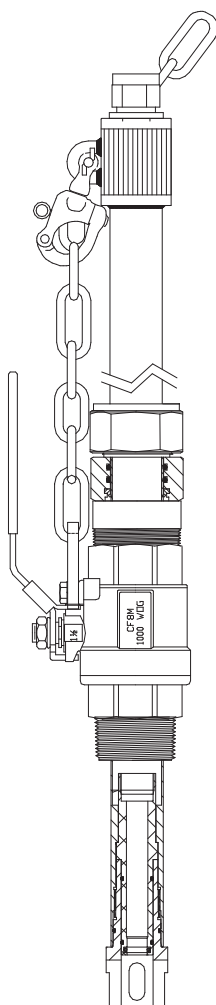
**InTrac® 787**

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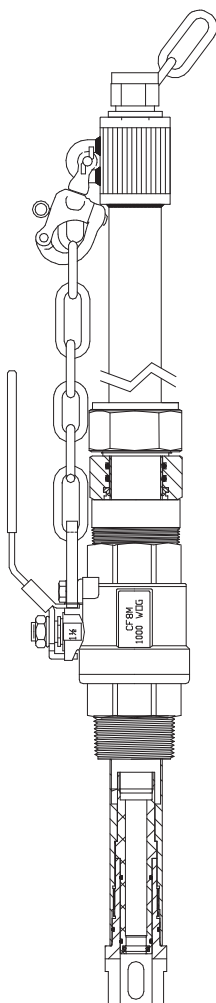
## Instruction manual

## Contents

<b>1</b>	<b>Introduction</b> .....	<b>5</b>
<b>2</b>	<b>Important remarks</b> .....	<b>6</b>
2.1	Checking the shipment .....	6
2.2	Notes on operating instructions .....	6
2.3	Normal use .....	7
2.4	Safety instructions .....	7
<b>3</b>	<b>Description of function</b> .....	<b>9</b>
<b>4</b>	<b>Installation/Startup and shutdown</b> .....	<b>10</b>
4.1	Preparation of the equipment.....	11
4.2	Fitting and installation works.....	11
4.2.1	Attaching the flushing lines .....	11
4.2.2	Fitting the electrode/sensor .....	12
4.4	Startup procedure.....	12
4.5	Shutdown procedure .....	12
<b>5</b>	<b>Maintenance</b> .....	<b>13</b>
5.1	Important information on maintenance .....	13
5.2	Replacing seals in contact with the process medium.....	13
<b>6</b>	<b>Trouble-shooting</b> .....	<b>14</b>
<b>7</b>	<b>Product specifications</b> .....	<b>15</b>
7.1	Standard equipment .....	15
7.2	Technical data .....	15
7.3	Spare parts list .....	16

# 1 Introduction

This operating instructions contains all the information needed for safe and proper use of the retractable housings **InTrac® 787**.

The operating instructions are intended for personnel entrusted with the operation and maintenance of the retractable housings. It is assumed that these persons are familiar with the equipment in which the retractable housing is installed.

Construction of the retractable housings employs leading-edge technology and complies with safety regulations currently in force (Mettler-Toledo GmbH as the manufacturer has been awarded **ISO 9001/EN29001** certification). Notwithstanding this, improper use could lead to hazards for the user or a third-party, and/or adverse effects on the plant or other equipment. Therefore, **the operating instructions must be read and understood** by the persons involved before work is started with the retractable housing.

In addition to these operating instructions, you should also be familiar with the following:

- all local safety regulations.
- all information and warnings in the documentation dealing with the products used together with the retractable housing (electrodes/sensors, control system, etc.).
- all safety regulations governing the equipment in which the retractable housing is installed.
- all information and warnings affixed to the retractable housing.

## 2 Important remarks

Please read through this section carefully. It contains several important points which will help you use the retractable housing in a safe and proper manner.

### 2.1 Checking the shipment

On receipt of the shipment, check immediately:

- the retractable housing and accessories for transport damage. Report any damage immediately to the carrier and your supplier.
- the type designation on the housing body.
- the shipment for completeness. Please notify your supplier immediately if the shipment is incomplete (see Section 7.1 "Standard equipment") or does not concur with the order and delivery documents.

### 2.2 Notes on operating instructions

#### Restriction

The information contained in this **operating instruction applies only to the InTrac® 787**. Detailed information on the electrodes and their installation is not given in these instructions. Such information can be found in the separate documentation accompanying these products. Where necessary, appropriate cross-references to such publications are given in these operating instructions.

#### Conventions

**This pictogram represents safety and hazard warnings** which, if ignored, could result in injuries to personnel and/or material damage.

#### Storage

Please keep these operating instructions in a safe place where they are always accessible. If these instructions are lost, please contact your supplier for a replacement.

## 2.3 Normal use

The InTrac® 787 is intended solely for measurement with the specified METTLER TOLEDO 12 mm pH and redox screwcap electrodes as well dissolved oxygen sensors.

Use the retractable housings solely for this purpose. Usage which differs from or exceeds this normal usage will be regarded as usage not for the intended purpose.

The manufacturer/supplier accepts no responsibility for any damage resulting from such improper usage. The risk is borne entirely by the user.

Normal use also includes:

- Compliance with the instructions, regulations and information in this operating instruction for the InTrac® 787 retractable housings.
- Adherence to the stipulated inspection and maintenance intervals.
- Correct maintenance of the retractable housing.
- Operation with due observance of the stipulated environmental and operating and the admissible mounting positions.

## 2.4 Safety instructions

The retractable housings InTrac® 787 should be installed, operated, maintained and, if need be, repaired only by personnel familiar with the housing and who are qualified for such work. The customer is responsible for ensuring that:

- only trained and authorized personnel work with the retractable housings.
- the operating instructions should be supplemented by in-house instructions regarding supervisory duties and obligation to report, work organization, personnel qualification, etc.
- the chain has to be always adjusted to the shortest possible length.

Ensure that the material specification of your retractable housing and of the seals satisfies the requirements of the process and is compatible with the (vessel) cleaning solutions used.

Before every startup, the retractable housing must be checked for:

- damage to the connections, fastening, etc.
- leaks

**A defective retractable housing must neither be installed nor put into service.** If it is suspected that hazard-free operation is no longer possible (e.g. housing damaged, connections leaking), the retractable housing and the equipment in which the retractable housing is installed must be shut down and secured against inadvertent operation.

Before dismantling the entire retractable housing or commencing any maintenance work on it, ensure that the equipment in which the housing is installed is in a safe condition (depressurize, empty, rinse, purge, vent, etc.).

**In the measuring position no maintenance work must be done.**

Only the maintenance and repair work described in these operating instructions may be performed on the retractable housings.

When changing faulty components, **use only original spare parts** obtainable from your supplier.

**No attachments or modifications** to the retractable housings and the accessories are allowed. The manufacturer/supplier accepts no responsibility for damage caused by unauthorized attachments and modifications or the incorporation of non-original spare parts. The risk is borne entirely by the user.

The rinse connections of the InTrac® 787 have a direct connection to the process when the ball valve is open. This means that you will have the effective process pressure on the rinsing tubes as well. Never do any manipulation on this connections while the ball valve is open.

### 3 Description of function

With the InTrac® 787 retractable housings, electrodes can be cleaned or replaced without interrupting the ongoing process. The process can be closed safely with the integrated ball valve. "The integrated flushing port allow a rinsing of the electrode in the housing.

The immersion tube is moved into the desired end position ("Measuring" or "Maintenance") by hand.

- In the "Measuring" position, the immersion tube is fully inserted in the process.
- In the "Maintenance" position, the electrode can be cleaned or removed/exchanged. The ball valve seals the housing against ingress of process medium.

## 4 Installation/Startup and shutdown

### 4.1 Preparation of the equipment

The housings are mounted and fixed on the vessel (reactor, tank, pipe, etc.) a flange connection or a 1" MNPT external (male) thread.

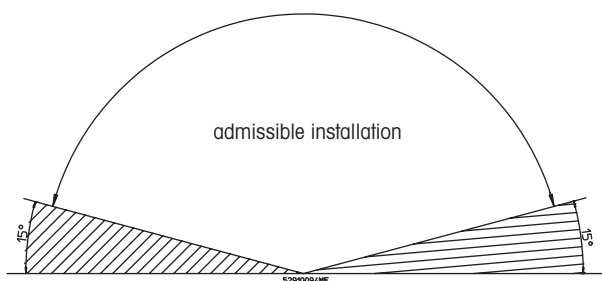
- The retractable housing should be mounted in a position which allows sufficient clearance for its correct functioning (correct measuring position in the medium) and maintenance (checks, installation and removal of the electrodes/sensors).
- Avoid mounting the retractable housing at exposed positions. If this is not possible, appropriate measures must be taken to protect against damage.

In order to guarantee correct function of the retractable housings, please pay attention to the following installation instructions:

- The retractable housing can be mounted vertically or in an inclined position.

**Caution!** In the case of inclined mounting, the angle of the housing must be equal to or greater than  $15^\circ$  above the horizontal.

### 4.2 Fitting and installation works

**Warning!** For all installation work described below, ensure that the equipment in which the housing is installed is in a non-hazardous condition (depressurized, empty, rinsed, purged, etc.).

#### 4.2.1 Attaching the flushing lines

##### Important information

- When the electrode/sensor is withdrawn, small amounts of the process medium adhere to the electrode/sensor tip and can enter the outflow line during the rinsing process. If the process medium is an environmentally harmful substance (toxic, corrosive, etc.), it is essential to comply with the local regulations governing the installation of wastewater services.
- Use only corrosion-resistant and chemically resistant materials for the rinsing water installation.

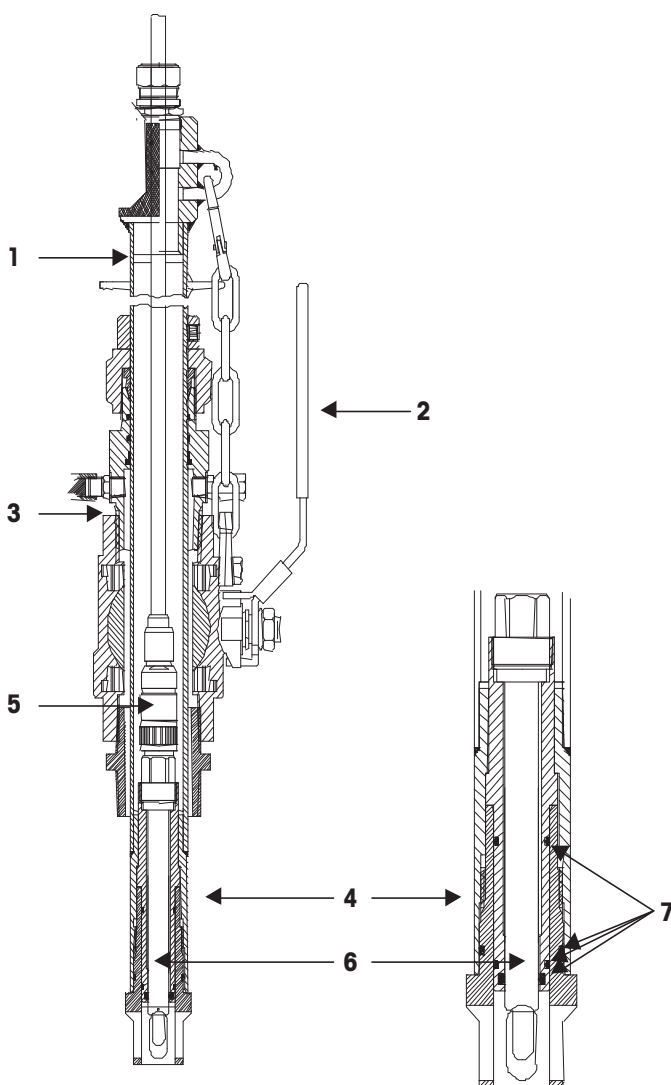
**Recommendation for proper operation and safe use**

Flushing lines must be equipped with valves which may only be opened when the retractable part of the housing is in the maintenance position. After the flushing and maintenance of the sensors, the flushing valves must be closed.

**Basic installation**

The basic installation is intended for cleaning/flushing of the electrodes/sensors when the housing is in the "Maintenance" position. In the optional accessories you will find Serto or Swagelok fitting to connect your flushing lines. The position of the Inlet and Outlet has no influence to the performance.

## 4.2.2 Fitting the electrode/sensor


1. Pull out the immersion tube.
2. Close the ball valve.
3. Disconnect the immersion tube right at the ball valve.
4. Disconnect the electrode holder by turning it counterclockwise.
5. Disconnect the electrode cable.
6. Remove the electrode.
7. Check washers and O-rings on electrodes for damage and replace if necessary.
8. Remove watering cap from electrode tip and rinse electrode/sensor with water.
9. Carefully insert the electrode into the electrode holder and screw in by hand as far as it will go.  
**Important:** use no tools!
10. Re-Install the complete housing.

### 4.3 Startup procedure

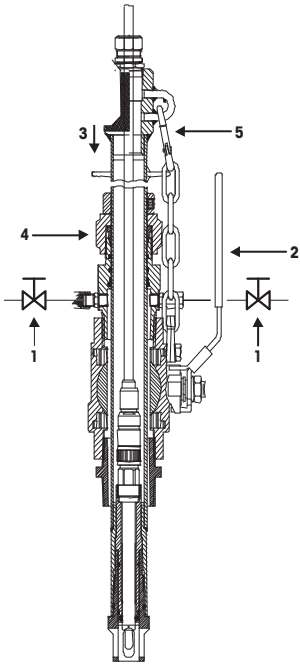
Before startup, all fitting and installation work must have been completed and the immersion tube must be secured to the ball valve with the safety chain!

1. **Warning!** Close the valves on both connections of the flushing chamber. Otherwise process media can escape while the ball valve is opened!

2. Open the ball valve.
3. Carefully push the immersion tube by hand into the process.
4. Secure immersion tube in this position by tightening the locking nut.
5. Shorten the safety chain to the shortest possible length.

### 4.4 Shutdown procedure

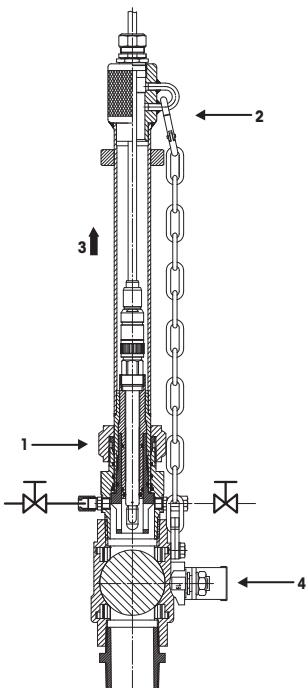1. Loosen the locking nut.
2. Hold immersion tube, then disconnect the safety chain.
3. Move immersion tube by hand to "Maintenance" position.
4. Close the ball valve.

5. **Caution!** Make sure that the process vessel or line is in a non-hazardous condition (depressurized, empty, rinsed, purged, etc.) when the housing has to be removed.

## 5 Maintenance

### 5.1 Important information on maintenance

**Attention!** Maintenance and service work on the housings may only be carried out by trained personnel.

**Warning!** It is possible that the process medium could harm the environment and your health (toxic, corrosive, etc.). Hence, ensure that the equipment is in a non-hazardous condition before you start with maintenance work.

- The electrode and housing must kept clean.
- Replace any damages seals or components of the housing
- Details in the maintenance of the electrode are contained in the electrode operation instructions.

### 5.2 Replacing seals in contact with the process medium

**All seals which come into contact with the process medium should always be replaced at least once a year for reasons of safety.**

With corrosive media, the seals may need to be changed at correspondingly shorter intervals.

**Warning!**

- Retractable housings may only be stripped down after having been completely dismantled.
- Check the quality of the seals and examine for possible damage. Ensure correct positioning.

## 6 Trouble-shooting

In this section you will find an overview of possible malfunctions which can occur during operation of the retractable housing, as well as information on how to rectify them.

<b>Malfunction</b>	<b>Cause</b>	<b>Remedial action</b>
Immersion tube cannot be inserted.	Ball valve is not completely opened.	Check the ball valve position.
Immersion tube cannot be pulled out.	Retention nut is not open.	Loosen the retention nut.
Immersion tube cannot be pulled out completely.	Ball valve is not completely opened.	Check the ball valve position.
Process liquid is flowing out of the rinsing chamber.	Rinsing chamber is not closed.	Close the outlet of the rinsing chamber with covers or with valves.
Process fluid is entering into the immersion tube.	O-ring defective.	Replace the O-ring in the electrode holder (order O-ring set)

## 7 Product specifications

### 7.1 Standard equipment

InTrac® 787 is supplied with:

- Immersion tube (100 mm or 300 mm)
- Electrode holder
- Ball valve
- Safety chain
- Electrode holder
- Threat connector 1 1/2"
- Nut to secure immersion tube

### 7.2 Technical data

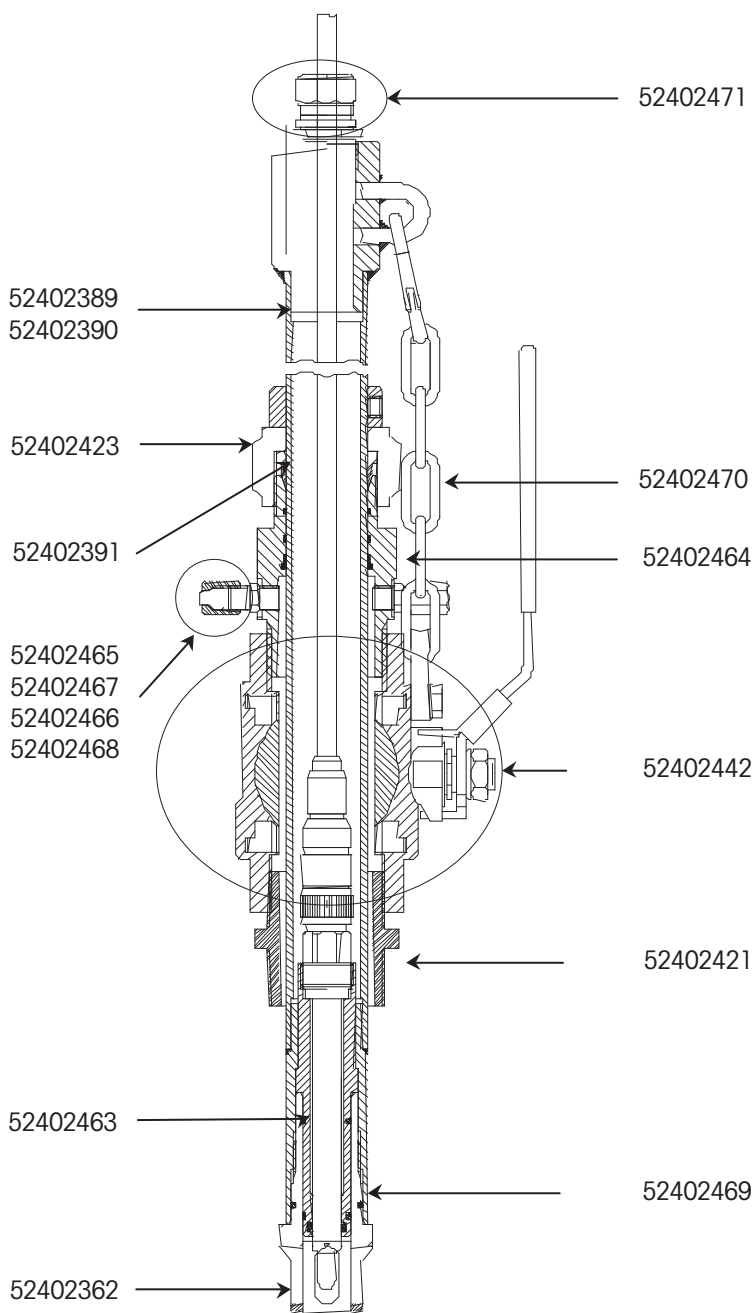
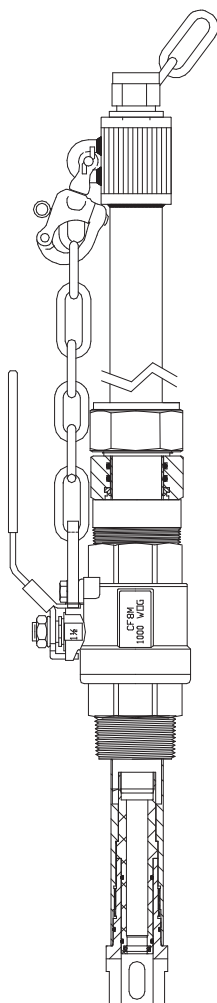
Dimensions	see dimensional drawing		
Maximum Pressure*	9 bar (130 psi)		
Maximum Temperature*	140°C ( 266 °F)		
Insertion Length	variable	0 to 100 (300) mm / 0" to 4(12)"	
Wetted Parts	Housing:	316L Stainless Steel (3.1B)	
	O-Rings:	Viton®	
	Ball valve:	Ball	316 (1.4401)
		Housing	CF-8M (1.4408)
	Sealing	PTFE	
Ball Valve	DN40, 1.5" NPT full port		

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- CH** Mettler-Toledo (Schweiz) AG, Im Langacher, Postfach, CH - 8606 Greifensee
- D** Mettler-Toledo GmbH, Prozeßanalytik, Ockerweg 3, D - 35396 Gießen
Analyse Industrielle S.A.S., 30, Boulevard Douaumont, BP 949, F - 75829 Paris Cedex 17
- CN** Mettler-Toledo Instruments (Shanghai) Co. Ltd.
- UK** Mettler-Toledo LTD, 64 Boston Road, Beaumont Leys, GB - Leicester LE4 1AW
- USA** Mettler-Toledo Ingold, Inc., 36 Middlesex Turnpike, Bedford, MA 01730, USA